

1 kb Plus DNA Ladder

Cat. no. 4992800/4992801

Storage: 4°C for 2 years (For long term storage, store at -20°C)

Concentration: About 0.1 mg DNA/ml

Product size

4992800	300 µl (50 lanes)
4992801	4 × 300 µl (200 lanes)

TIANGEN BIOTECH (BEIJING) CO., LTD.

[HTTP://WWW.TIANGEN.COM/EN](http://www.tiangen.com/en)

The product is used for research only, neither intended for the diagnosis, or treatment of a disease, nor for the food, or cosmetics etc.

Description

The 1 kb Plus DNA Ladder consists of twelve double-stranded DNA bands that are suitable as molecular weight standards for agarose gel electrophoresis.

Since this product is premixed with 1× Loading Buffer, it is convenient to load 3~6 µl of this product to agarose gel directly.

The ladder is composed of twelve individual DNA fragments (in base pairs): 10000, 8000, 5000, 2000, 1600, 1000, 7000, 500, 400, 300, 200 and 100. It contains one intense band (1000 bp) for easy orientation. To load 6 µl of the product will provide 50 ng per DNA fragment with one intense fragment of 100 ng.

Buffer Contents

10 mM Tris-HCl (pH 8.4)
 10 mM EDTA
 0.02% Bromophenol Blue
 5% Glycerol


Protocol

1. Directly load 3~6 µl of this product to the agarose gel (load 1 µl product per 1 mm gel lane. The loading volume can be increased appropriately for wide lanes).
2. Electrophoresis condition: 1.0% agarose gel, 4~10 v/cm voltage.
3. DNA marker can be visualized using ethidium bromide under UV light.

Recommendations

1. Directly use, do not require heating.
2. For ideal results, please use new agarose gel and change electrophoresis buffer in time.

1 kb Plus DNA Ladder

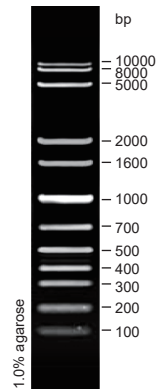

6 µl product,
 10 cm gel length,
 6 v/cm voltage,
 0.5 × TBE Buffer